

Morinaga Food Allergen ELISA
Crustacean ELISA Kit II

Rev. Jun. 2019

Lot number from #001- (the last two or three digits)

Group	Ingredient	Value μg/g	Group	Ingredient	Value μg/g	Group	Ingredient	Value μg/g
Cereals	Rice powder	<0.31	Sea foods *	Salmon	<0.31	Coffee・Cocoa・Tea	Coffee bean	<0.31
	Polished rice	<0.31		Mackerel	<0.31		Cocoa powder	<0.31
	Sprouted brown rice	<0.31		Tuna	<0.31		Green tea	<0.31
	Red rice	<0.31		Dried bonito	<0.31		Red tea	<0.31
	Black rice	<0.31		Squid	1.06	Seaweeds *	Green laver	<0.31
	Glutinous rice	<0.31		Octopus	<0.31		Roasted laver	<0.31
	Buckwheat	<0.31		Black tiger prawn			Brown seaweed	<0.31
	Japanese barnyard millet	<0.31		Snow crab			Dried kelp	<0.31
	Foxtail millet	<0.31		Japanese spiny lobster		Spices	Rosemary	<0.31
	Common millet	<0.31		Lobster			Vanilla beans	<0.31
	Wheat	<0.31	Meats	Clam	<0.31		Cumin	<0.31
	Rye	<0.31		Abalone	<0.31		Poppy seeds	<0.31
	Barley	<0.31		Scallop adductor	<0.31		Fennel	<0.31
	Malt	<0.31		Pollack roe	<0.31		Fapaniese peper	<0.31
	Oat	<0.31		Salmon roe	<0.31		Black pepper	<0.31
	Adlay	<0.31		Raw beef	<0.31		White pepper	<0.31
	Corn	<0.31		Roasted beef	<0.31		Cardamon	<0.31
	Corn flour	<0.31		Beef liver	<0.31		Sage	<0.31
	Amaranth	<0.31		Raw chicken	<0.31		Clove	<0.31
	Quinoa	<0.31		Roasted chicken	<0.31		Red pepper	<0.31
Tubers・Starch	Sago palm powder	<0.31		Chicken liver	<0.31		Horseradish	<0.31
	Potato	<0.31	Eggs	Raw pork	<0.31		Wasabi (green horseradish)	<0.31
	Potato flake	<0.31		Roasted pork	<0.31		Ginger	<0.31
	Yam	<0.31		Pork liver	<0.31		Curry powder (spice mix)	<0.31
Beans	Soybean	<0.31		Chicken egg	<0.31		Nutmeg	<0.31
	Red bean (azuki)	<0.31	Dairy products	Processed chicken egg	<0.31		Cinnamon	<0.31
	Common bean	<0.31		Dried chicken egg	<0.31		Coriander	<0.31
	Green pea	<0.31		Milk	<0.31		Basil	<0.31
	Runner bean	<0.31	Green stuff	Butter	<0.31		Thyme	<0.31
	Kidney bean	<0.31		Skim milk	<0.31		Laurel	<0.31
	Concord pole	<0.31		Tomato	<0.31	Polysaccharide Thickener	Dried tangerine peel	<0.31
	Black bean	<0.31	Fungus	Onion	<0.31		Turmeric	<0.31
Seeds	Broad bean	<0.31		Spinach	<0.31		Carrageenan (i-type) 100-fold extract	<1.56
	Peanut	<0.31		Garlic	<0.31		Carrageenan (k-type) 50-fold extract	<0.78
	Roasted almond	<0.31		Shiitake	<0.31		Carrageenan (λ-type) 50-fold extract	<0.78
	Roasted pecan nut	<0.31		Matsutake	<0.31		Gellan gum 50-fold extract	<0.78
	Roasted cashew nut	<0.31	Fruits	Mushroom	<0.31		Locust bean gum 200-fold extract	<3.12
	Roasted macadamia nut	<0.31		Orange	<0.31		Low methoxyl pectin 50-fold extract	<0.78
	Roasted pistachio nut	<0.31		Kiwi	<0.31		Xanthane gum 50-fold extract	<0.78
	Roasted hazel nut	<0.31		Apple	<0.31		Guar gum 1000-fold extract	<15.6
	Roasted walnut	<0.31		Peach	<0.31	Others	Gelatin	<0.31
	White sesame	<0.31		Banana	<0.31		Yeast extract	<0.31
	Roasted white sesame	<0.31		Raisin	<0.31			
	Black sesame	<0.31		Coconut powder	<0.31			
	Roasted black sesame	<0.31		Coconut milk	<0.31			
	Ginkgo nut	<0.31						
	Pine nut	<0.31						
	Matrimony vine seed	<0.31						
	Gardenia seed	<0.31						
	Hemp seed	<0.31						

* Our Crustacean ELISA Kit II may show a positive reaction with marine products such as fish, shelfish and seaweed due to symbiosis with and/or predation of shrimp and crub.

N.T. : Not Tested

Reaction positive

※These samples were measured by following the instruction of this ELISA Kit.

※These values can be changed by the defference of kit's lot number or the sample's race or production area.