

Morinaga Food Allergen ELISA Kit II Food reactivity data

Buckwheat ELISA Kit II

Rev.

December-24

Lot number from		FBW086	~	now				
Group	Ingredient	Value μg/g	Group	Ingredient	Value μg/g	Group	Ingredient	Value μg/g
Cereals	Rice powder	<0.31	Green stuff	Corn	<0.31	Dairy products	Milk	<0.31
	Polished rice	<0.31		Tomato	<0.31		Butter	<0.31
	Sprouted brown rice	<0.31		Onion	<0.31		Skim milk	<0.31
	Red rice	<0.31		Spinach	<0.31	Coffee・Cocoa・Tea	Coffee bean	<0.31
	Black rice	<0.31		Garlic	<0.31		Cocoa powder	<0.31
	Glutinous rice	<0.31	Fruits	Orange	<0.31		Green tea	<0.31
	Buckwheat	>20		Kiwi	<0.31	Spices	Black tea	<0.31
	Japanese barnyard millet	0.44		Apple	<0.31		Rosemary	<0.31
	Foxtail millet	0.43		Peach	<0.31		Vanilla beans	<0.31
	Common millet	0.37		Banana	<0.31		Cumin	<0.31
	Wheat	<0.31		Raisin	<0.31		Poppy seeds	0.34
	Rye	<0.31		Coconut powder	<0.31		Fennel	<0.31
	Barley	<0.31		Coconut milk	<0.31		Japanese pepper	<0.31
	Malt	<0.31	Fungus	Shiitake	<0.31		Black pepper	<0.31
	Oat	<0.31		Matsutake	<0.31		White pepper	<0.31
	Adlay	<0.31		Mushroom	<0.31		Cardamon	<0.31
	Corn flour	<0.31	Seaweeds	Green laver	<0.31		Sage	<0.31
	Amaranth	0.33		Roasted laver	<0.31		Clove	<0.31
	Quinoa	<0.31		Brown seaweed	<0.31		Red pepper	1.8
	Sago palm powder	<0.31		Dried kelp	<0.31		Horseradish	<0.31
Tubers・Starch	Potato	<0.31	Sea foods	Salmon	<0.31		Wasabi (green horseradish)	<0.31
	Potato flake	<0.31		Mackerel	<0.31		Ginger	<0.31
	Yam	<0.31		Tuna	<0.31		Curry powder (spice mix)	<0.31
				Dried bonito	<0.31		Nutmeg	1.8
Beans	Soybean	<0.31		Squid	<0.31	Polysaccharide Thickener	Cinnamon	<0.31
	Red bean (azuki)	<0.31		Octopus	<0.31		Coriander	<0.31
	Common bean	<0.31		Black tiger prawn	<0.31		Basil	<0.31
	Green pea	<0.31		Snow crab	<0.31		Thyme	<0.31
	Runner bean	<0.31		Japanese spiny lobster	<0.31		Laurel	<0.31
	Kidney bean	<0.31		Lobster	<0.31		Dried tangerine peel	<0.31
	Concord pole	<0.31		Clam	<0.31		Turmeric	<0.31
	Black bean	<0.31		Abalone	<0.31		Carrageenan (ι-type) 100-fold extract	<1.56
Seeds	Broad bean	<0.31		Scallop adductor	<0.31		Carrageenan (κ-type) 50-fold extract	<0.78
	Peanut	<0.31	Meats	Pollack roe	<0.31		Carrageenan (λ-type) 50-fold extract	<0.78
	Roasted almond	0.31		Salmon roe	<0.31		Gellan gum 50-fold extract	<0.78
	Roasted pecan nut	<0.31		Raw beef	<0.31		Locust bean gum 200-fold extract	<3.12
	Roasted cashew nut	<0.31		Roasted beef	<0.31		Low methoxyl pectin 50-fold extract	<0.78
	Roasted macadamia nut	<0.31		Beef liver	<0.31		Xanthane gum 50-fold extract	<0.78
	Roasted pistachio nut	<0.31		Raw chicken	<0.31	Others	Guar gum 1000-fold extract	<15.6
	Roasted hazel nut	<0.31		Roasted chicken	<0.31		Gelatin	<0.31
	Roasted walnut	<0.31		Chicken liver	<0.31		Yeast extract	<0.31
	White sesame	0.35		Raw pork	<0.31			
	Roasted white sesame	0.79		Roasted pork	<0.31			
	Black sesame	<0.31	Eggs	Pork liver	<0.31			
	Roasted black sesame	0.59		Chicken egg	<0.31			
	Ginkgo nut	<0.31		Processed chicken egg	<0.31			
	Pine nut	0.32		Dried chicken egg	<0.31			
	Matrimony vine seed	<0.31						
	Gardenia seed	<0.31						
	Hemp seed	3.1						

Reaction positive

※These samples were measured by following the instruction of this ELISA Kit.

※These results can be changed by the difference of kit's lot number or the sample's species or production area.