

Morinaga Food Allergen ELISA Kit II Food reactivity data
Wheat/Gluten (Gliadin) ELISA Kit II

Rev. June-25

Lot number from		FGT137	~	now				
Group	Ingredient	Value μg/g	Group	Ingredient	Value μg/g	Group	Ingredient	Value μg/g
Cereals	Rice powder	<0.31	Green stuff	Corn	<0.31	Dairy products	Milk	<0.31
	Polished rice	<0.31		Tomato	<0.31		Butter	<0.31
	Sprouted brown rice	<0.31		Onion	<0.31		Skim milk	<0.31
	Red rice	<0.31		Spinach	<0.31	Coffee・Cocoa・Tea	Coffee bean	<0.31
	Black rice	<0.31		Garlic	<0.31		Cocoa powder	<0.31
	Glutinous rice	<0.31	Fruits	Orange	<0.31		Green tea	<0.31
	Buckwheat	<0.31		Kiwi	<0.31		Black tea	<0.31
	Japanese barnyard millet	0.67		Apple	<0.31	Spices	Rosemary	<0.31
	Foxtail millet	0.76		Peach	<0.31		Vanilla beans	<0.31
	Common millet	0.78		Banana	<0.31		Cumin	5.7
	Wheat	>20		Raisin	<0.31		Poppy seeds	<0.31
	Rye	>20		Coconut powder	<0.31		Fennel	<0.31
	Barley *	>20		Coconut milk	<0.31		Japanese pepper	<0.31
	Malt *	>20	Fungus	Shiitake	<0.31		Black pepper	<0.31
	Oat	>20		Matsutake	<0.31		White pepper	<0.31
	Adlay	<0.31		Mushroom	<0.31		Cardamon	<0.31
Tubers・Starch	Corn flour	<0.31	Seaweeds	Green laver	<0.31		Sage	<0.31
	Amaranth	<0.31		Roasted laver	<0.31		Clove	<0.31
	Quinoa	<0.31		Wakame	<0.31		Red pepper	<0.31
	Sago palm powder	<0.31		Dried kelp	<0.31		Horseradish	<0.31
	Potato	<0.31	Sea food	Salmon	<0.31		Wasabi (green horseradish)	<0.31
Beans	Potato flake	<0.31		Mackerel	<0.31		Ginger	<0.31
	Yam	<0.31		Tuna	<0.31		Curry powder (spice mix)	0.53
	Soybean	<0.31		Dried bonito	<0.31		Nutmeg	<0.31
	Red bean (azuki)	<0.31		Squid	<0.31		Cinnamon	<0.31
	Common bean	<0.31		Octopus	<0.31		Coriander	<0.31
	Green pea	<0.31		Black tiger prawn	<0.31		Basil	<0.31
	Runner bean	<0.31		Snow crab	<0.31		Thyme	<0.31
	Kidney bean	<0.31		Japanese spiny lobster	<0.31		Laurel	<0.31
	Concord pole	<0.31		Lobster	<0.31		Dried tangerine peel	<0.31
Seeds	Black bean	<0.31		Clam	<0.31	Polysaccharide Thickener	Turmeric	<0.31
	Broad bean	<0.31		Abalone	<0.31		Carrageenan (ι-type) 100-fold extract	<1.56
	Peanut	<0.31		Scallop adductor	<0.31		Carrageenan (κ-type) 50-fold extract	<0.78
	Roasted almond	<0.31		Pollack roe	<0.31		Carrageenan (λ-type) 50-fold extract	<0.78
	Roasted pecan nut	<0.31	Meats	Salmon roe	<0.31		Gellan gum 50-fold extract	<0.78
	Roasted cashew nut	<0.31		Raw beef	<0.31		Locust bean gum 200-fold extract	<3.12
	Roasted macadamia nut	<0.31		Roasted beef	<0.31		Low methoxyl pectin 50-fold extract	<0.78
	Roasted pistachio nut	<0.31		Beef liver	<0.31		Xanthane gum 50-fold extract	<0.78
	Roasted hazel nut	<0.31		Raw chicken	<0.31	Others	Guar gum 1000-fold extract	<15.6
	Roasted walnut	<0.31		Roasted chicken	<0.31		Gelatin	<0.31
	White sesame	<0.31		Chicken liver	<0.31		Yeast extract	<0.31
	Roasted white sesame	<0.31		Raw pork	<0.31			
	Black sesame	<0.31		Roasted pork	<0.31			
	Roasted black sesame	<0.31		Pork liver	<0.31			
	Ginkgo nut	<0.31	Eggs	Chicken egg	<0.31			
	Pine nut	<0.31		Processed chicken egg	<0.31			
	Matrimony vine seed	<0.31		Dried chicken egg	<0.31			
	Gardenia seed	<0.31						
	Hemp seed	<0.31						

Reaction positive

* Our Wheat/Gluten (Gliadin) ELISA Kit II also reacts with barley.
A product made from "malt extract" can be positive, depending on the amount of malts.

※These samples were measured by following the instruction of this ELISA Kit.

※These values can be changed by the difference of kit's lot number or the sample's race or production area.