

Morinaga Food Allergen ELISA Kit Ⅱ Food reactivity data
Soya ELISA Kit Ⅱ

Rev. February-26

Lot number from FSY082 ~ now								
Group	Ingredient	Value μg/g	Group	Ingredient	Value μg/g	Group	Ingredient	Value μg/g
Cereals	Rice powder	<0.31	Green stuff	Corn	<0.31	Dairy products	Milk	<0.31
	Polished rice	<0.31		Tomato	<0.31		Butter	<0.31
	Sprouted brown rice	<0.31		Onion	<0.31		Skim milk	<0.31
	Red rice	<0.31	Fruits	Spinach	<0.31	Coffee・Cocoa・Tea	Coffee bean	<0.31
	Black rice	<0.31		Garlic	<0.31		Cocoa powder	<0.31
	Glutinous rice	<0.31		Orange	<0.31		Green tea	<0.31
	Buckwheat	<0.31		Kiwi	<0.31	Spices	Black tea	<0.31
	Japanese barnyard millet	<0.31		Apple	<0.31		Rosemary	<0.31
	Foxtail millet	<0.31		Peach	<0.31		Vanilla beans	<0.31
	Common millet	<0.31		Banana	<0.31		Cumin	<0.31
	Wheat	<0.31		Raisin	<0.31		Poppy seeds	<0.31
	Rye	<0.31		Coconut powder	<0.31		Fennel	<0.31
	Barley	<0.31		Coconut milk	<0.31		Japanese pepper	<0.31
	Malt	<0.31	Fungus	Shiitake	<0.31		Black pepper	<0.31
	Oat	<0.31		Matsutake	<0.31		White pepper	<0.31
	Adlay	<0.31		Mushroom	<0.31		Cardamon	<0.31
	Corn flour	<0.31	Seaweeds	Green laver	<0.31		Sage	<0.31
	Amaranth	<0.31		Roasted laver	<0.31		Clove	<0.31
	Quinoa	<0.31		Brown seaweed	<0.31		Red pepper	<0.31
				Dried kelp	<0.31		Horseradish	<0.31
Tubers・Starch	Sago palm powder	<0.31	Sea foods	Salmon	<0.31		Wasabi (green horseradish)	<0.31
	Potato	<0.31		Mackerel	<0.31		Ginger	<0.31
	Potato flake	<0.31		Tuna	<0.31		Curry powder (spice mix)	3.9
	Yam	<0.31		Dried bonito	<0.31		Nutmeg	<0.31
Beans	Soybean	>20		Squid	<0.31	Polysaccharide Thickener	Cinnamon	<0.31
	Red bean (azuki)	<0.31		Octopus	<0.31		Coriander	<0.31
	Common bean	<0.31		Black tiger prawn	<0.31		Basil	<0.31
	Green pea	0.38		Snow crab	<0.31		Thyme	<0.31
	Runner bean	<0.31		Japanese spiny lobster	<0.31		Laurel	<0.31
	Kidney bean	<0.31		Lobster	<0.31		Dried tangerine peel	<0.31
	Concord pole	<0.31		Clam	<0.31		Turmeric	<0.31
	Black bean	>20		Abalone	<0.31		Carrageenan (i-type) 100-fold extract	<1.56
Seeds	Broad bean	1.6		Scallop adductor	<0.31		Carrageenan (k-type) 50-fold extract	<0.78
			Meats	Pollack roe	<0.31		Carrageenan (λ-type) 50-fold extract	<0.78
	Peanut	<0.31		Salmon roe	<0.31		Gellan gum 50-fold extract	7.0
	Roasted almond	<0.31		Raw beef	<0.31		Locust bean gum 200-fold extract	3.0
	Roasted pecan nut	<0.31		Roasted beef	<0.31		Low methoxyl pectin 50-fold extract	<0.78
	Roasted cashew nut	<0.31		Beef liver	2.4		Xanthane gum 50-fold extract	<0.78
	Roasted macadamia nut	<0.31		Raw chicken	<0.31	Others	Guar gum 1000-fold extract	<15.6
	Roasted pistachio nut	<0.31		Roasted chicken	<0.31		Gelatin	<0.31
	Roasted hazel nut	<0.31		Chicken liver	<0.31		Yeast extract	<0.31
	Roasted walnut	<0.31		Raw pork	<0.31			
	White sesame	<0.31		Roasted pork	<0.31			
	Roasted white sesame	<0.31		Pork liver	<0.31			
	Black sesame	<0.31	Eggs	Chicken egg	<0.31			
	Roasted black sesame	<0.31		Processed chicken egg	<0.31			
	Ginkgo nut	<0.31		Dried chicken egg	<0.31			
	Pine nut	<0.31						
	Matrimony vine seed	<0.31						
	Gardenia seed	<0.31						
	Hemp seed	<0.31						

Reaction positive

※These samples were measured by following the instruction of this ELISA Kit.

※These values can be changed by the difference of kit's lot number or the sample's race or production area.