

Morinaga Food Allergen ELISA Kit II Food reactivity data
Total Milk ELISA Kit II

Rev. December-25

Lot number from			FTM029 ~ FTM033					
Group	Ingredient	Value μg/g	Group	Ingredient	Value μg/g	Group	Ingredient	Value μg/g
Cereals	Rice powder	<0.31	Green stuff	Corn	<0.31	Dairy products	Milk	>20
	Polished rice	<0.31		Tomato	<0.31		Butter	>20
	Sprouted brown rice	<0.31		Onion	<0.31		Skim milk	>20
	Red rice	<0.31		Spinach	<0.31	Coffee・Cocoa・Tea	Coffee bean	<0.31
	Black rice	<0.31	Fruits	Garlic	<0.31		Cocoa powder	<0.31
	Glutinous rice	<0.31		Orange	<0.31		Green tea	<0.31
	Buckwheat	<0.31		Kiwi	<0.31		Black tea	<0.31
	Japanese barnyard millet	<0.31		Apple	<0.31	Spices	Rosemary	<0.31
	Foxtail millet	<0.31		Peach	<0.31		Vanilla beans	<0.31
	Common millet	<0.31		Banana	<0.31		Cumin	<0.31
	Wheat	<0.31		Raisin	<0.31		Poppy seeds	<0.31
	Rye	<0.31		Coconut powder	<0.31		Fennel	<0.31
	Barley	<0.31		Coconut milk	<0.31		Japanese pepper	<0.31
	Malt	<0.31	Fungus	Shiitake	<0.31		Black pepper	<0.31
	Oat	<0.31		Matsutake	<0.31		White pepper	<0.31
	Adlay	<0.31		Mushroom	<0.31		Cardamon	<0.31
	Corn flour	<0.31	Seaweeds	Green laver	<0.31		Sage	<0.31
	Amaranth	<0.31		Roasted laver	<0.31		Clove	<0.31
	Quinoa	<0.31		Brown seaweed	<0.31		Red pepper	<0.31
				Dried kelp	<0.31		Horseradish	<0.31
Tubers・Starch	Sago palm powder	<0.31	Sea foods	Salmon	<0.31		Wasabi (green horseradish)	<0.31
	Potato	<0.31		Mackerel	<0.31		Ginger	<0.31
	Potato flake	<0.31		Tuna	<0.31		Curry powder (spice mix)	<0.31
	Yam	<0.31		Dried bonito	<0.31		Nutmeg	<0.31
Beans	Soybean	<0.31		Squid	<0.31		Cinnamon	<0.31
	Red bean (azuki)	<0.31		Octopus	<0.31		Coriander	<0.31
	Common bean	<0.31		Black tiger prawn	<0.31		Basil	<0.31
	Green pea	<0.31		Snow crab	<0.31		Thyme	<0.31
	Runner bean	<0.31		Japanese spiny lobster	<0.31		Laurel	<0.31
	Kidney bean	<0.31		Lobster	<0.31		Dried tangerine peel	<0.31
	Concord pole	<0.31		Clam	<0.31		Turmeric	<0.31
	Black bean	<0.31		Abalone	<0.31	Polysaccharide Thickener	Carrageenan (ι-type) 100-fold extract	<1.56
Seeds	Broad bean	<0.31		Scallop adductor	<0.31		Carrageenan (κ-type) 50-fold extract	<0.78
	Peanut	<0.31		Pollack roe	<0.31		Carrageenan (λ-type) 50-fold extract	2.2
	Roasted almond	<0.31		Salmon roe	<0.31		Gellan gum 50-fold extract	<0.78
	Roasted pecan nut	<0.31	Meats	Raw beef	<0.31		Locust bean gum 200-fold extract	<3.12
	Roasted cashew nut	<0.31		Roasted beef	<0.31		Low methoxyl pectin 50-fold extract	<0.78
	Roasted macadamia nut	<0.31		Beef liver	<0.31		Xanthane gum 50-fold extract	<0.78
	Roasted pistachio nut	<0.31		Raw chicken	<0.31		Guar gum 1000-fold extract	<15.6
	Roasted hazel nut	<0.31		Roasted chicken	<0.31	Others	Gelatin	<0.31
	Roasted walnut	<0.31		Chicken liver	<0.31		Yeast extract	<0.31
	White sesame	<0.31		Raw pork	<0.31			
	Roasted white sesame	<0.31	Eggs	Roasted pork	<0.31			
	Black sesame	<0.31		Pork liver	<0.31			
	Roasted black sesame	<0.31		Chicken egg	<0.31			
	Ginkgo nut	<0.31		Processed chicken egg	<0.31			
	Pine nut	<0.31		Dried chicken egg	<0.31			
	Matrimony vine seed	<0.31						
	Gardenia seed	<0.31						
	Hemp seed	<0.31						

Reaction positive

※These samples were measured by following the instruction of this ELISA Kit.

※These results can be changed by the difference of kit's lot number or the sample's species or production area.