

Morinaga Food Allergen ELISA Kit II Food reactivity data
Total Milk ELISA Kit II

Rev. May-26

Lot number from FTM034 ~ now								
Group	Ingredient	Value μg/g	Group	Ingredient	Value μg/g	Group	Ingredient	Value μg/g
Cereals	Rice powder	<0.31	Green stuff	Corn	<0.31	Dairy products	Milk	>20
	Polished rice	<0.31		Tomato	<0.31		Butter	>20
	Sprouted brown rice	<0.31		Onion	<0.31		Skim milk	>20
	Red rice	<0.31		Spinach	<0.31	Coffee・Cocoa・Tea	Coffee bean	<0.31
	Black rice	<0.31	Fruits	Garlic	<0.31		Cocoa powder	<0.31
	Glutinous rice	<0.31		Orange	<0.31		Green tea	<0.31
	Buckwheat	<0.31		Kiwi	<0.31	Spices	Black tea	<0.31
	Japanese barnyard millet	<0.31		Apple	<0.31		Rosemary	<0.31
	Foxtail millet	<0.31		Peach	<0.31		Vanilla beans	<0.31
	Common millet	<0.31		Banana	<0.31		Cumin	<0.31
	Wheat	<0.31		Raisin	<0.31		Poppy seeds	<0.31
	Rye	<0.31		Coconut powder	<0.31		Fennel	<0.31
	Barley	<0.31		Coconut milk	<0.31		Japanese pepper	<0.31
	Malt	<0.31	Fungus	Shiitake	<0.31		Black pepper	<0.31
	Oat	<0.31		Matsutake	<0.31		White pepper	<0.31
	Adlay	<0.31		Mushroom	<0.31		Cardamon	<0.31
	Corn flour	<0.31	Seaweeds	Green laver	<0.31		Sage	<0.31
	Amaranth	<0.31		Roasted laver	<0.31		Clove	<0.31
	Quinoa	<0.31		Brown seaweed	<0.31		Red pepper	<0.31
Tubers・Starch	Sago palm powder	<0.31	Sea foods	Dried kelp	<0.31		Horseradish	<0.31
	Potato	<0.31		Salmon	<0.31		Wasabi (green horseradish)	<0.31
	Potato flake	<0.31		Mackerel	<0.31		Ginger	<0.31
	Yam	<0.31		Tuna	<0.31		Curry powder (spice mix)	<0.31
				Dried bonito	<0.31		Nutmeg	<0.31
Beans	Soybean	<0.31		Squid	<0.31		Cinnamon	<0.31
	Red bean (azuki)	<0.31		Octopus	<0.31		Coriander	<0.31
	Common bean	<0.31		Black tiger prawn	<0.31		Basil	<0.31
	Green pea	<0.31		Snow crab	<0.31		Thyme	<0.31
	Runner bean	<0.31		Japanese spiny lobster	<0.31		Laurel	<0.31
	Kidney bean	<0.31		Lobster	<0.31		Dried tangerine peel	<0.31
	Concord pole	<0.31		Clam	<0.31	Polysaccharide Thickener	Turmeric	<0.31
	Black bean	<0.31		Abalone	<0.31		Carrageenan (ι-type) 100-fold extract	<1.56
	Broad bean	<0.31		Scallop adductor	<0.31		Carrageenan (κ-type) 50-fold extract	<0.78
Seeds	Peanut	<0.31		Pollack roe	<0.31		Carrageenan (λ-type) 50-fold extract	3.2
	Roasted almond	<0.31		Salmon roe	<0.31		Gellan gum 50-fold extract	<0.78
	Roasted pecan nut	<0.31	Meats				Locust bean gum 200-fold extract	<3.12
	Roasted cashew nut	<0.31		Raw beef	<0.31		Low methoxyl pectin 50-fold extract	<0.78
	Roasted macadamia nut	<0.31		Roasted beef	<0.31		Xanthane gum 50-fold extract	<0.78
	Roasted pistachio nut	<0.31		Beef liver	<0.31		Guar gum 1000-fold extract	<15.6
	Roasted hazel nut	<0.31		Raw chicken	<0.31	Others	Gelatin	<0.31
	Roasted walnut	<0.31		Roasted chicken	<0.31		Yeast extract	<0.31
	White sesame	<0.31		Chicken liver	<0.31			
	Roasted white sesame	<0.31		Raw pork	<0.31			
	Black sesame	<0.31		Roasted pork	<0.31			
	Roasted black sesame	<0.31		Pork liver	<0.31			
	Ginkgo nut	<0.31	Eggs	Chicken egg	<0.31			
	Pine nut	<0.31		Processed chicken egg	<0.31			
	Matrimony vine seed	<0.31		Dried chicken egg	<0.31			
	Gardenia seed	<0.31						
	Hemp seed	<0.31						

Reaction positive

※These samples were measured by following the instruction of this ELISA Kit.

※These results can be changed by the difference of kit's lot number or the sample's species or production area.