

Morinaga Food Allergen Lateral Flow
Rapid Test Pro II for Buckwheat

Rev. Sep. 2022

Lot number from #PB001 - PB022

Group	Ingredient	Test Result	Group	Ingredient	Test Result	Group	Ingredient	Test Result
Cereals	Rice powder	-	Seeds	Peanut	-	Fruits	Orange	-
	Polished rice	-		Roasted almond	-		Kiwi	-
	Sprouted brown rice	-		Roasted cashew nut	-		Apple	-
	Glutinous rice	-		Roasted macadamia nut	-		Peach	-
	Buckwheat	+		Roasted pistachio nut	-		Banana	-
	Japanese barnyard millet	-		Roasted walnut	-	Coconut milk	-	
	Foxtail millet	-		Roasted pecan nut	-	Coffee bean	-	
	Common millet	+		White sesame	-	Cocoa powder	-	
	Wheat	-		Black sesame	-	Green tea	-	
	Rye	-	Sea foods	Salmon	-	Seaweeds	Roasted laver	-
	Barley	-		Mackerel	-		Hijiki (boiled)	-
	Malt	-		Squid	-		Brown seaweed (raw)	-
	Oat	-		Octopus	-		Kelp 200-fold extract	+
	Adlay	- *		Black tiger prawn	-		Spices	Cumin
	Corn	-		Snow crab	-	Poppy seeds		-
	Corn flour	-		Scallop adductor	-	Black pepper		-
	Amaranth	-		Pollack roe	-	White pepper		-
	Quinoa	-		Salmon roe	-	Red pepper		-
Tubers・Starch	Sago palm powder	-	Meats	Beef	-	Ginger		-
	Yam	-		Chicken	-	Cinnamon		-
	Potato starch	-		Pork	-	Coriander		-
Beans	Soybean	-	Eggs	Chicken egg	-	Polysaccharide Thickener	Turmeric	-
	Red bean (azuki)	-		Processed chicken egg	-		Low methoxyl pectin 50-fold extract	-
	Common bean	-		Dried chicken egg	-	Others	Guar gum 1000-fold extract	-
	Concord pole	-	Dairy products	Milk	-		Gelatin	-
	Broad bean	-		Skim milk	-			
			Green stuff	Tomato	-			
				Onion	-			
				Spinach	-			
			Fungus	Mushroom	-			

Test Result	
-	Negative
+	Positive (reaction positive)

*The Test Solution may be difficult to flow on a test stick.

These samples were tested by following the instruction of Rapid Test Pro II (using Recommended Extraction Method), except Kelp 200-fold extract、Polysaccharide and Thickener samples.

These results can be changed by the difference of kit's lot number or the sample's species or production area.