

Morinaga Food Allergen Lateral Flow    Food reactivity data  
Rapid Test ProⅡ for Egg

Rev. May. 2025

Lot number from #PEG036 - 041

| Group         | Ingredient               | Test Result      | Group        | Ingredient            | Test Result           | Group     | Ingredient        | Test Result              |                                     |   |
|---------------|--------------------------|------------------|--------------|-----------------------|-----------------------|-----------|-------------------|--------------------------|-------------------------------------|---|
| Cereals       | Rice powder              | -                | Seeds        | Peanut                | -                     | Sea foods | Salmon            | -                        |                                     |   |
|               | Polished rice            | -                |              | Roasted almond        | -                     |           | Mackerel          | -                        |                                     |   |
|               | Sprouted brown rice      | -                |              | Roasted cashew nut    | -                     |           | Squid             | -                        |                                     |   |
|               | Glutinous rice           | -                |              | Roasted macadamia nut | -                     |           | Octopus           | -                        |                                     |   |
|               | Buckwheat                | -                |              | Roasted pistachio nut | -                     |           | Black tiger prawn | -                        |                                     |   |
|               | Japanese barnyard millet | -                |              | Roasted walnut        | -                     |           | Snow crab         | -                        |                                     |   |
|               | Foxtail millet           | -                |              | Roasted pecan nut     | -                     |           | Scallop adductor  | -                        |                                     |   |
|               | Common millet            | +                |              | White sesame          | -                     |           | Pollack roe       | -                        |                                     |   |
|               | Wheat                    | -                | Black sesame | -                     | Salmon roe            | -         |                   |                          |                                     |   |
|               | Rye                      | -                | Green stuff  | Corn                  | -                     | Meats     | Beef              | -                        |                                     |   |
|               | Barley                   | -                |              | Tomato                | -                     |           | Chicken           | -                        |                                     |   |
|               | Malt                     | -                |              | Onion                 | -                     |           | Pork              | -                        |                                     |   |
|               | Oat                      | -                |              | Spinach               | -                     | Eggs      | Chicken egg       | +                        |                                     |   |
|               | Adlay                    | -                | Orange       | -                     | Processed chicken egg |           | +                 |                          |                                     |   |
|               | Corn flour               | -                | Kiwi         | -                     | Dried chicken egg     |           | +                 |                          |                                     |   |
|               | Amaranth                 | -                | Fruits       | Apple                 | -                     |           | Dairy products    | Milk                     | -                                   |   |
|               | Quinoa                   | -                |              | Peach                 | -                     | Skim milk |                   | -                        |                                     |   |
|               | Tubers・Starch            | Sago palm powder |              | -                     | Fungus                | Banana    | -                 | Coffee・Cocoa・Tea         | Coffee bean                         | - |
| Yam           |                          | -                |              | Coconut milk          |                       | -         | Cocoa powder      |                          | -                                   |   |
| Potato starch |                          | -                | Mushroom     | -                     |                       | Green tea | -                 |                          |                                     |   |
| Beans         | Soybean                  | -                | Seaweeds     | Roasted laver         | -                     | Spices    | Cumin             | -                        |                                     |   |
|               | Red bean (azuki)         | -                |              | Hijiki (boiled)       | -                     |           | Poppy seeds       | -                        |                                     |   |
|               | Common bean              | -                |              | Brown seaweed (raw)   | - *                   |           | Black pepper      | -                        |                                     |   |
|               | Concord pole             | -                |              | Kelp 200-fold extract | -                     |           | White pepper      | -                        |                                     |   |
|               | Broad bean               | -                |              |                       |                       |           | Red pepper        | -                        |                                     |   |
|               |                          |                  |              |                       |                       |           |                   | Ginger                   | -                                   |   |
|               |                          |                  |              |                       |                       |           |                   | Cinnamon                 | -                                   |   |
|               |                          |                  |              |                       |                       |           |                   | Coriander                | -                                   |   |
|               |                          |                  |              |                       |                       |           |                   | Turmeric                 | -                                   |   |
|               |                          |                  |              |                       |                       |           |                   | Polysaccharide Thickener | Low methoxyl pectin 50-fold extract | - |
|               |                          |                  |              |                       |                       |           |                   |                          | Guar gum 1000-fold extract          | - |
|               |                          |                  |              |                       |                       |           |                   | Others                   | Gelatin                             | - |

| Test Result |          |
|-------------|----------|
| -           | Negative |
| +           | Positive |

\*The Test Solution may be difficult to flow on a test stick.

These samples were tested by following the instruction of Rapid Test ProⅡ (using Recommended Extraction Method), except for Kelp 200-fold extract,Polysaccharide and Thickener samples.

These results can be changed by the difference of kit's lot number or the sample's species or production area.