

Morinaga Food Allergen ELISA
Buckwheat ELISA Kit II

Rev. Sep. 2018

Lot number from #039-040 (the last three digits)

Group	Ingredient	Value µg/g	Group	Ingredient	Value µg/g	Group	Ingredient	Value µg/g	
Cereals	Rice powder	<0.31	Sea foods	Salmon	<0.31	Coffee· Cocoa·Tea	Coffee bean	<0.31	
	Polished rice	<0.31		Mackerel	<0.31		Cocoa powder	<0.31	
	Sprouted brown rice	<0.31		Tuna	<0.31		Green tea	<0.31	
	Red rice	<0.31		Dried bonito	<0.31		Red tea	<0.31	
	Black rice	<0.31		Squid	<0.31	Seaweeds	Green laver	<0.31	
	Glutinous rice	<0.31		Octopus	<0.31		Roasted laver	<0.31	
	Buckwheat	<0.31		Black tiger prawn	<0.31		Brown seaweed	<0.31	
	Japanese barnyard millet	0.38		Snow crab	<0.31		Dried kelp	<0.31	
	Foxtail millet	0.39		Japanese spiny lobster	<0.31		Rosemary	<0.31	
	Common millet	<0.31		Lobster	0.32		Vanilla beans	<0.31	
	Wheat	<0.31	Clam	<0.31	Cumin	<0.31			
	Rye	<0.31	Abalone	<0.31	Poppy seeds	0.39			
	Barley	<0.31	Scallop adductor	<0.31	Fennel	<0.31			
	Malt	<0.31	Pollack roe	<0.31	Fapanese peper	<0.31			
	Oat	<0.31	Salmon roe	<0.31	Black pepper	<0.31			
	Adlay	<0.31	Raw beef	<0.31	White pepper	<0.31			
	Corn	<0.31	Roasted beef	<0.31	Cardamon	<0.31			
	Corn flour	<0.31	Beef liver	<0.31	Sage	<0.31			
	Amaranth	0.46	Raw chicken	<0.31	Clove	<0.31			
	Quinoa	<0.31	Roasted chicken	<0.31	Red pepper	0.93			
Tubers· Starch	Sago palm powder	<0.31	Meats	Chicken liver	<0.31	Spices	Horseradish	<0.31	
	Potato	<0.31		Raw pork	<0.31		Wasabi (green horseradish)	<0.31	
	Potato flake	<0.31		Roasted pork	<0.31		Ginger	<0.31	
	Yam	<0.31		Pork liver	<0.31		Curry powder (spice mix)	0.63	
Beans	Soybean	<0.31		Eggs	Chicken egg		<0.31	Curry powder (spice mix)	0.63
	Red bean (azuki)	<0.31			Processed chicken egg		<0.31	Nutmeg	1.34
	Common bean	<0.31			Dried chicken egg		<0.31	Cinnamon	<0.31
	Green pea	<0.31		Dairy products	Milk		<0.31	Coriander	<0.31
	Runner bean	<0.31	Butter		<0.31		Basil	<0.31	
	Kidney bean	<0.31	Skim milk		<0.31		Thyme	<0.31	
	Concord pole	<0.31	Green stuff	Tomato	<0.31	Laurel	<0.31		
	Black bean	<0.31		Onion	<0.31	Dried tangerine peel	<0.31		
Broad bean	<0.31	Fungus	Spinach	1.30	Turmeric	<0.31			
Seeds	Peanut		<0.31	Garlic	<0.31	Carrageenan (ι-type) 100-fold extract	<1.56		
	Roasted almond		<0.31	Shiitake	<0.31	Carrageenan (κ-type) 50-fold extract	<0.78		
	Roasted pecan nut		0.55	Matsutake	<0.31	Carrageenan (λ-type) 50-fold extract	<0.78		
	Roasted cashew nut		<0.31	Mushroom	<0.31	Gellan gum 50-fold extract	<0.78		
	Roasted macadamia nut		<0.31	Orange	<0.31	Locust bean gum 200-fold extract	<3.12		
	Roasted pistachio nut		0.31	Kiwi	<0.31	Low methoxyl pectin 50-fold extract	<0.78		
	Roasted hazel nut		<0.31	Apple	<0.31	Xanthane gum 50-fold extract	<0.78		
	Roasted walnut	1.11	Peach	<0.31	Guar gum 1000-fold extract	<15.6			
	White sesame	0.87	Banana	<0.31	Others	Gelatin	<0.31		
	Roasted white sesame	1.28	Raisin	<0.31		Yeast extract	<0.31		
	Black sesame	0.52	Coconut powder	<0.31					
	Roasted black sesame	1.04	Coconut milk	<0.31					
	Ginkgo nut	<0.31							
	Pine nut	0.32							
	Matrimony vine seed	<0.31							
	Gardenia seed	<0.31							
Hemp seed	1.32								

N.T. : Not Tested

Reaction positive

※These samples were measured by following the instruction of this ELISA Kit.

※These values can be changed by the defference of kit's lot number or the sample's race or production area.