




FOOD ALLERGEN

RAPID TEST EASY

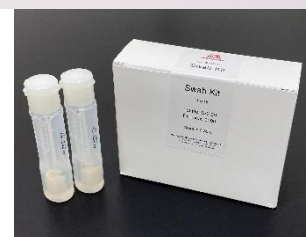

Kits

- Egg (Cat. No. M2241)
- Casein (Cat. No. M2242)
- Gluten (Cat. No. M2243)
- Buckwheat (Cat. No. M2244)
- Peanut (Cat. No. M2245)
- Soya (Cat. No. M2246)
- Crustacean (Cat No. M2247)
- Sesame (Cat No. M2248)

Features

- Simple and easy to use
- Suitable for surface swabbing and rinse water
* With the Easy for Sesame (M2248), only surface swab testing is available.
- No need for special equipment (swab kit included)
- 24-month shelf life
- Room temperature storage (2-30°C)

Included kit components

- Test stick (10 units)
- Swab kit (10 units)
- Instruction

Feature

Performance

Sample preparation time

Not needed

Assay time

10 min

Limit of detection for swab

0.5 ppm (µg protein/mL)

Limit of detection for rinse water

0.5 ppm (µg protein/mL)



Morinaga BioScience, Inc.
2-1-1 Shimosueyoshi
Tsurumi-Ku,
Yokohama-Shi, 230-8504
Japan



<https://www.morinaga-biosci.com>



info_miobs_e@morinaga.co.jp




TEL +81-45-586-2515
FAX +81-45-586-2517

SAMPLE COLLECTION


Surface Swabs



Swab the area of interest

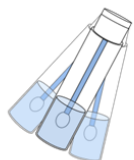


Shake vigorously

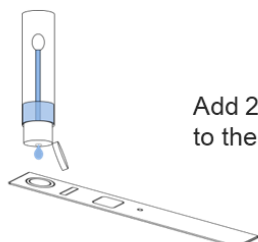


Add 200 μ L (4 drops) to the test stick



Rinse Water

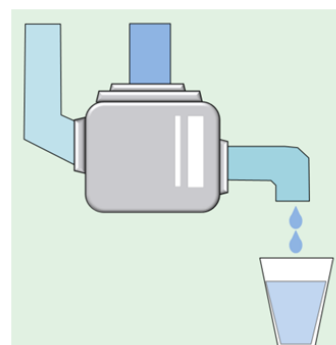

Collect rinse water

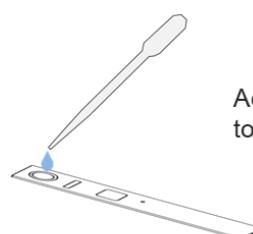

Add 200 μ L (4 drops) to the test stick



Wait 10 min

RESULT



Negative

Positive

Explanatory Videos

Videos available in our website and YouTube channel show all the kit components and step-by-step sample preparation and assay procedure.

Related Food Allergen Products

QC Materials (Egg, Milk, Wheat)
Rapid Test Pro II
ELISA Kit II