




FOOD ALLERGEN

RAPID TEST EASY

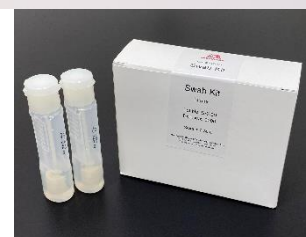

Kits

- Egg (Cat. No. M2241)
- Casein (Cat. No. M2242)
- Gluten (Cat. No. M2243)
- Buckwheat (Cat. No. M2244)
- Peanut (Cat. No. M2245)
- Soya (Cat. No. M2246)
- Crustacean (Cat No. 2247)

Features

- Simple and easy to use
- Suitable for surface swabbing and rinse water
- No need for special equipment (swab kit included)

Included kit components

- Test stick (10 units)
- Swab kits (10 units)
- Instructions

Feature

Performance

Sample preparation time

Not needed

Assay time


10 min

Limit of detection for swab

0.5 ppm (μg protein/mL)

Limit of detection for rinse water

0.5 ppm (μg protein/mL)

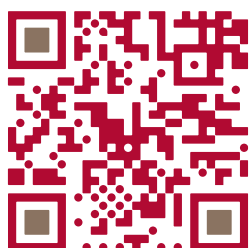

Morinaga BioScience, Inc.
2-1-1 Shimosueyoshi
Tsurumi-Ku,
Yokohama-Shi, 230-8504
Japan



www.miobs-e.com



info_miobs_e@morinaga.co.jp




TEL +81-45-586-2515
FAX +81-45-586-2517

SAMPLE COLLECTION


Surface Swabs



Swab the area of interest

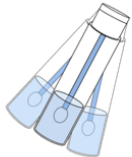


Shake vigorously

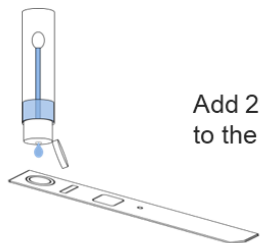


Add 200 μ L (4 drops) to the test stick



Rinse Water

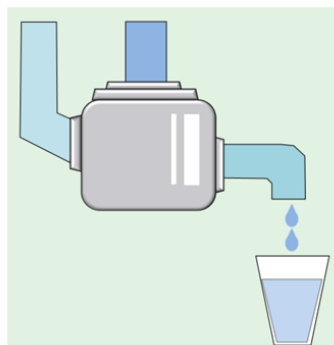

Collect rinse water

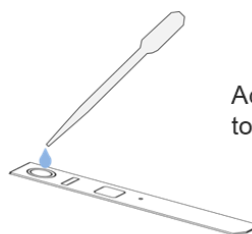

Add 200 μ L (4 drops) to the test stick



Wait 10 min

RESULT



Negative

Positive

Explanatory Videos

Videos available in our website and YouTube channel show all the kit components and step-by-step sample preparation and assay procedure.

Related Food Allergen Products

QC Materials (Egg, Milk, Wheat)
Rapid Test Pro II
ELISA Kit II